



Closed Loop and Ceramic Nozzles

BD Aria's


Lift the
Aria cover to
view ceramic
nozzle.

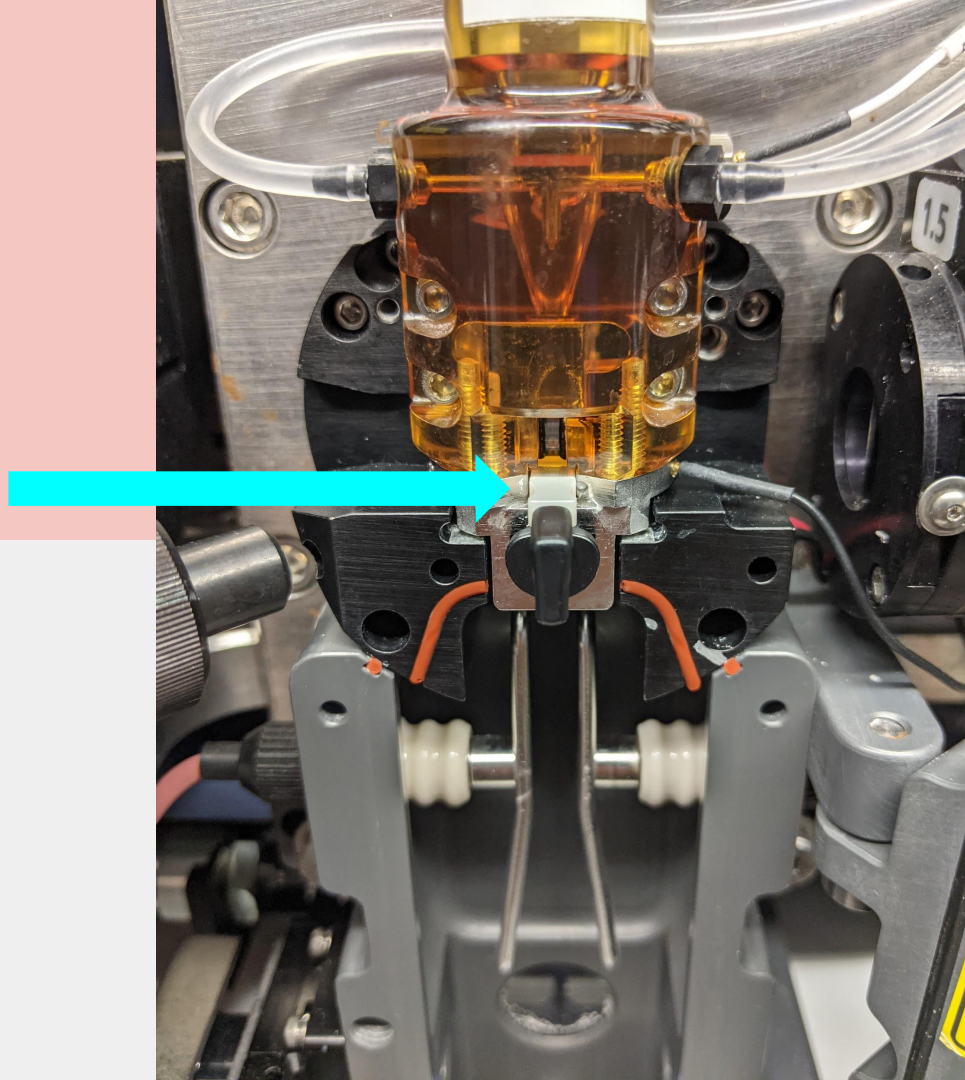


To remove the ceramic nozzle, turn the black knob counter clockwise to 9 o'clock.

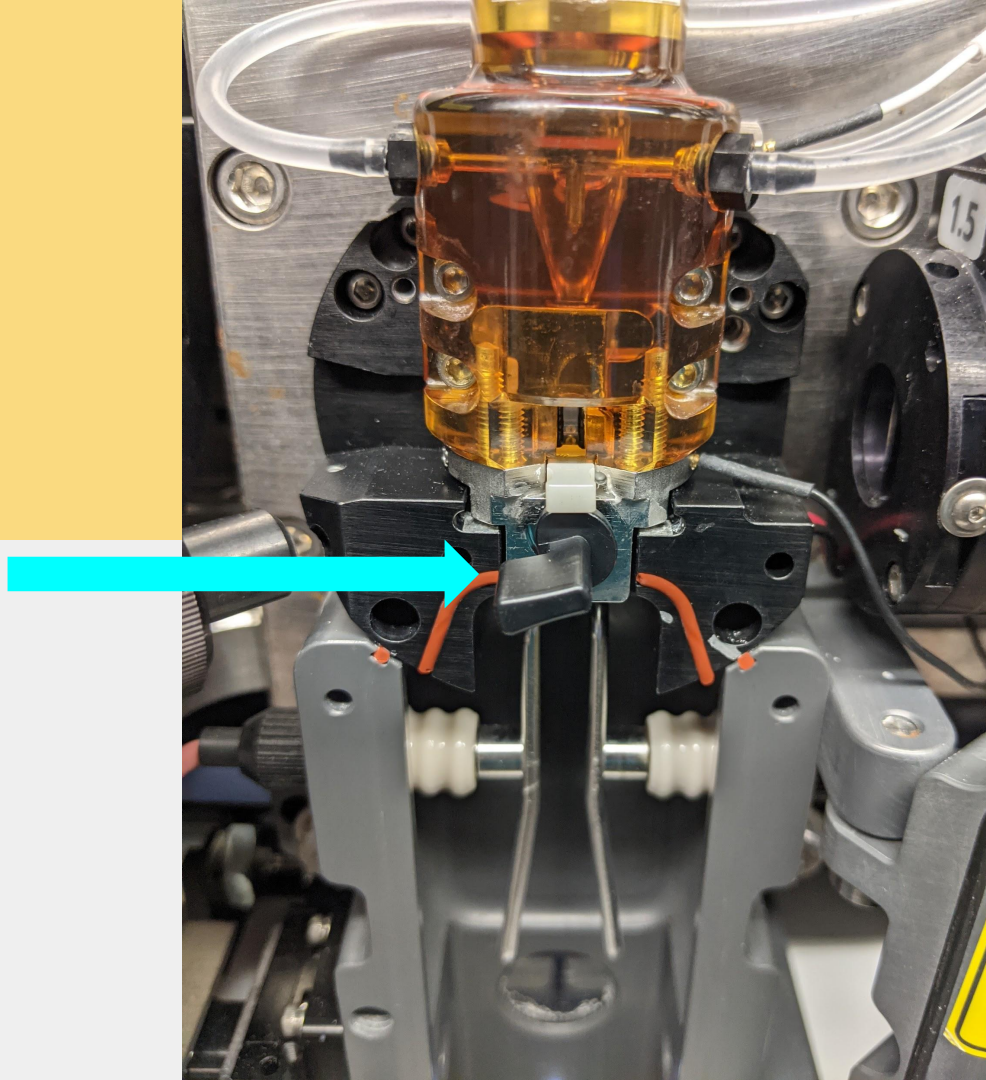


Pull the ceramic nozzle straight out.

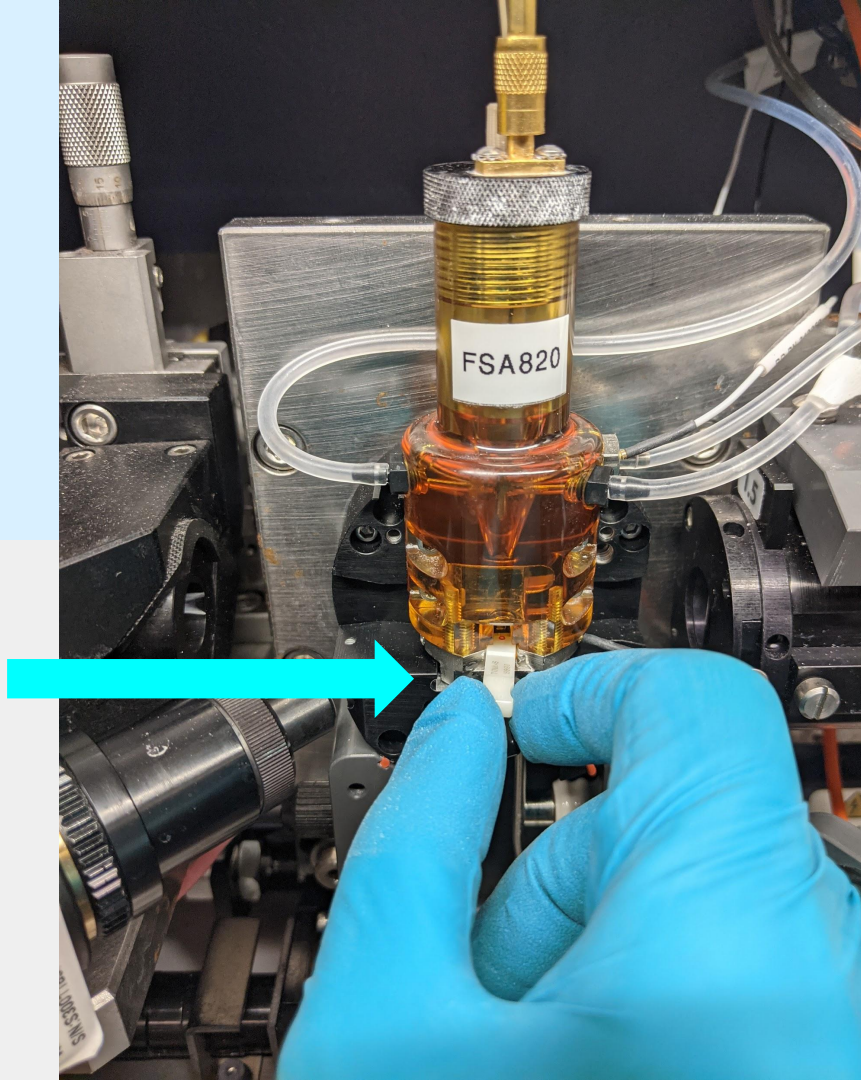


No nozzle in the
Quartz Cuvette.

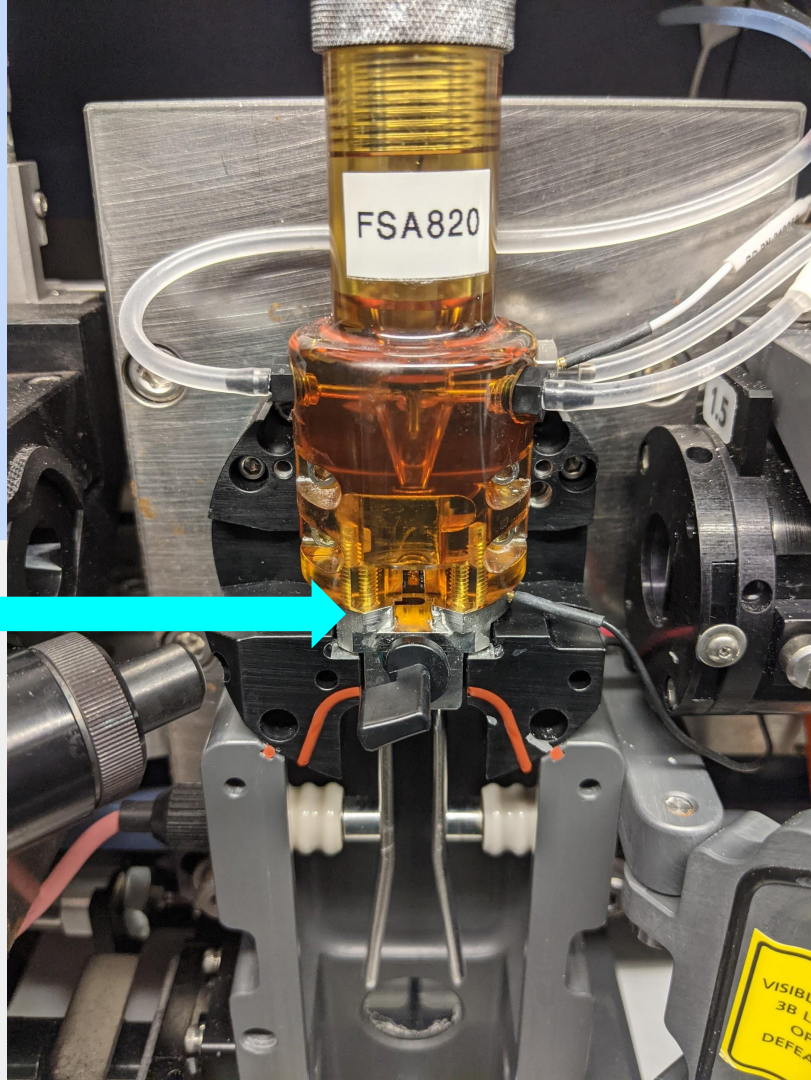



Put the ceramic nozzle in the 15ml water tube marked NOZZLE and put it in the sort chamber.

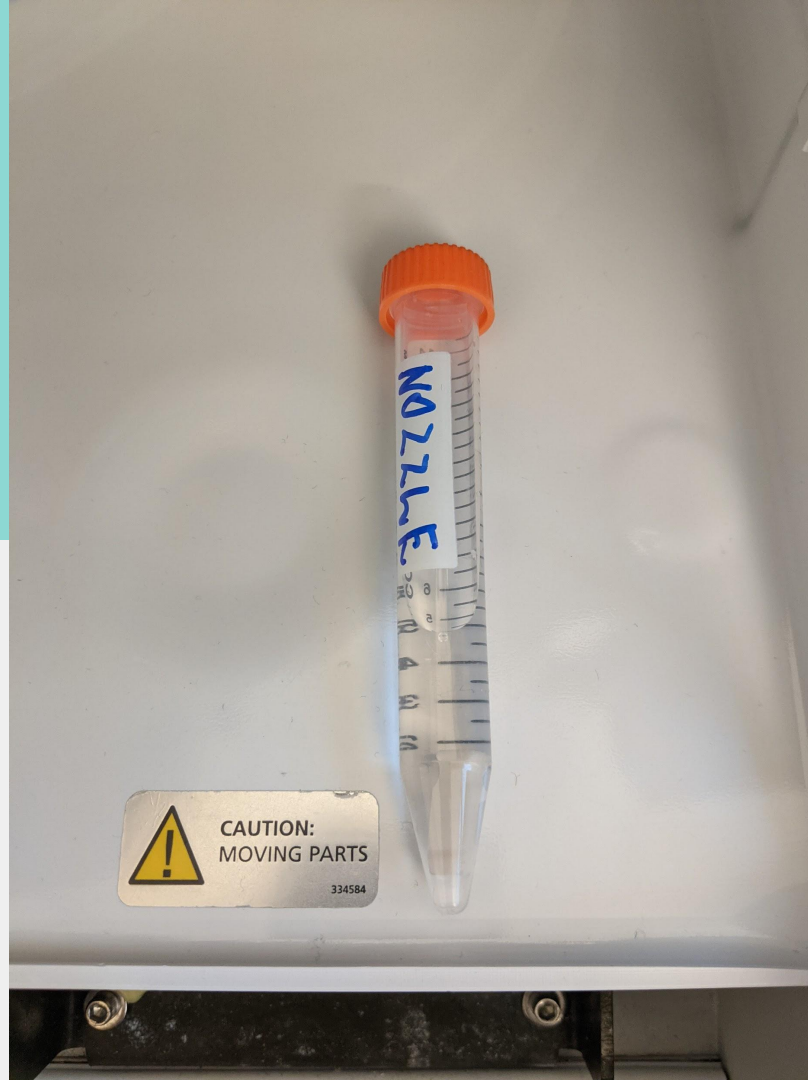


Locate the Closed Loop Nozzle to the right of the Quartz Cuvette. It has the same shape as the ceramic nozzle with a brown nut and vacuum tubing attached to it.

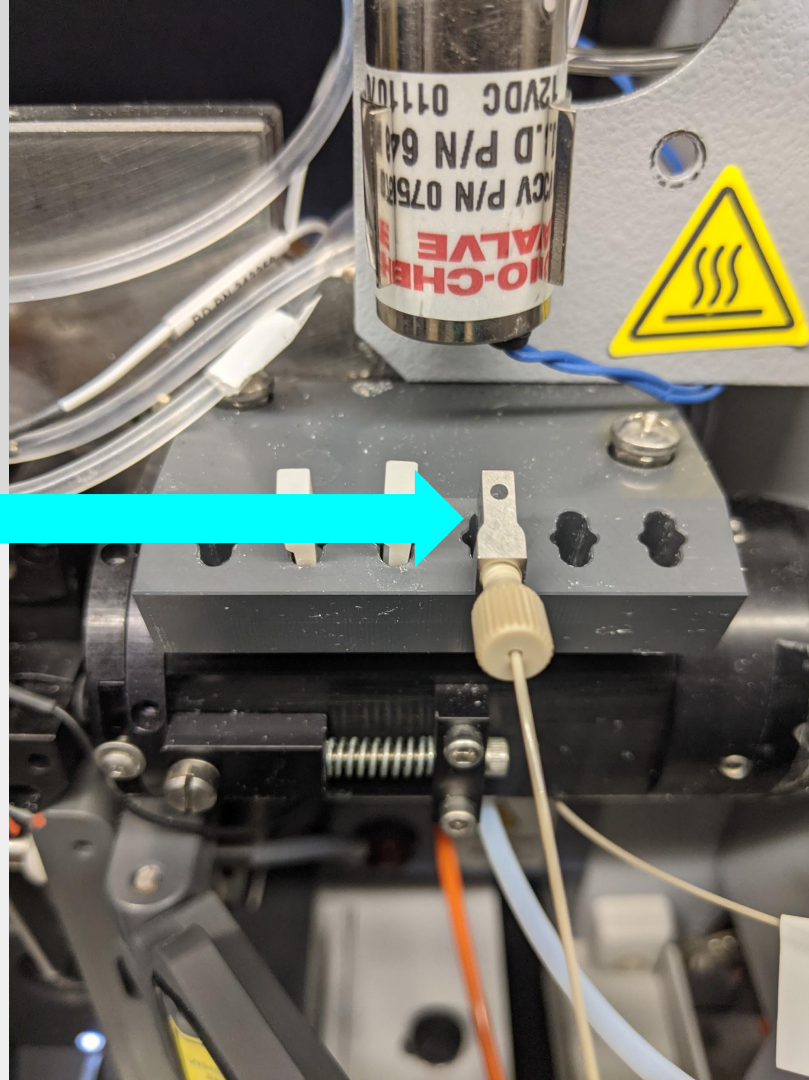


Notice the
Closed Loop
Nozzle is labeled
as such.

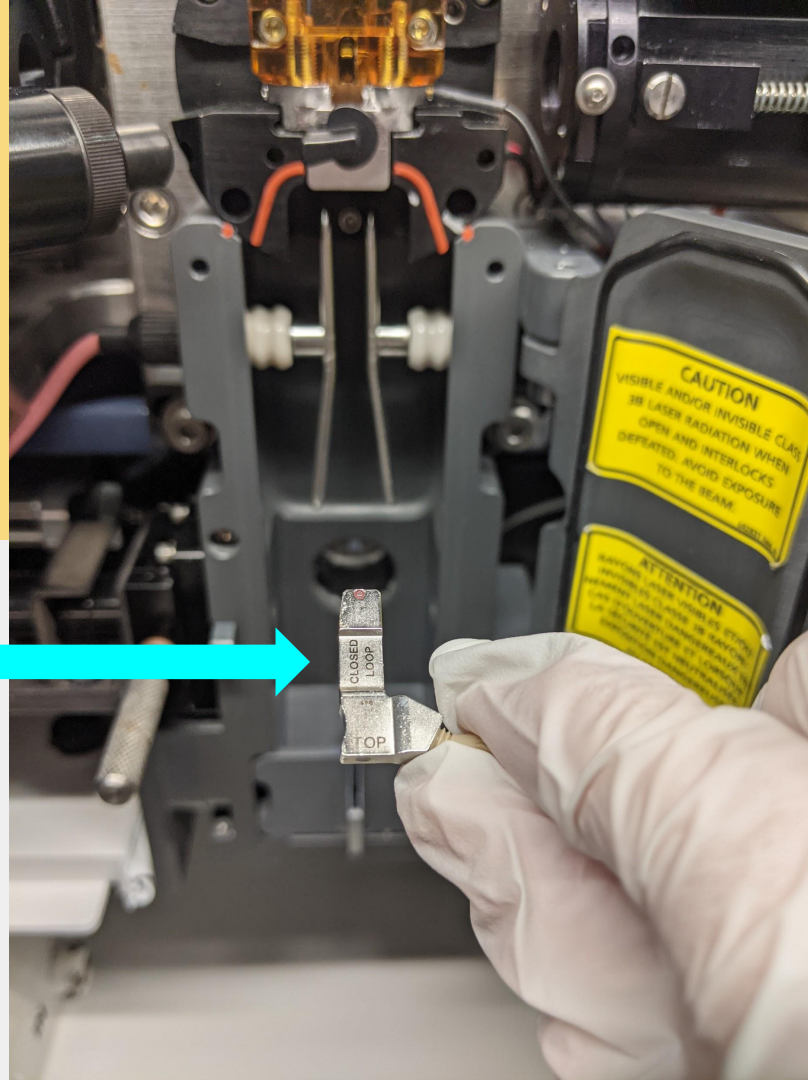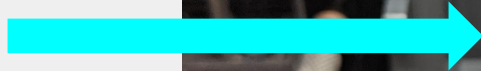

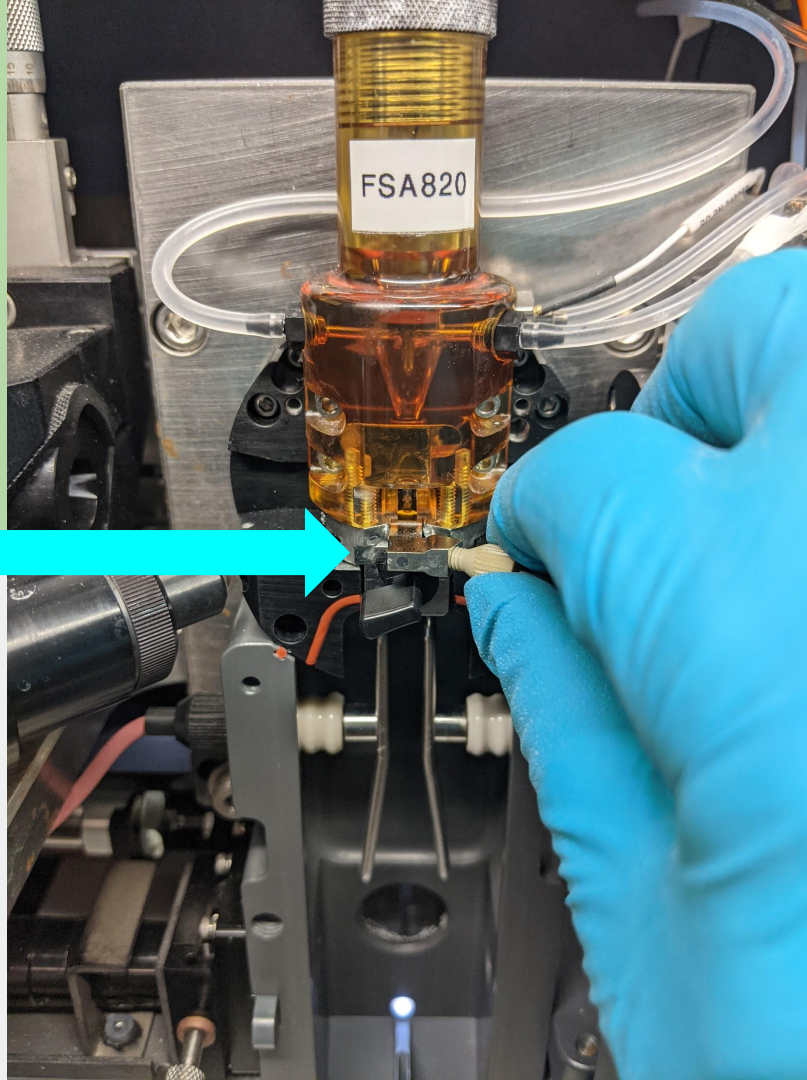

The Closed Loop Nozzle red o ring and label faces upwards when inserted into the Quartz Cuvette.

Turn the black knob clockwise to 12 o'clock to seat the red o ring up against the Quart Cuvette.

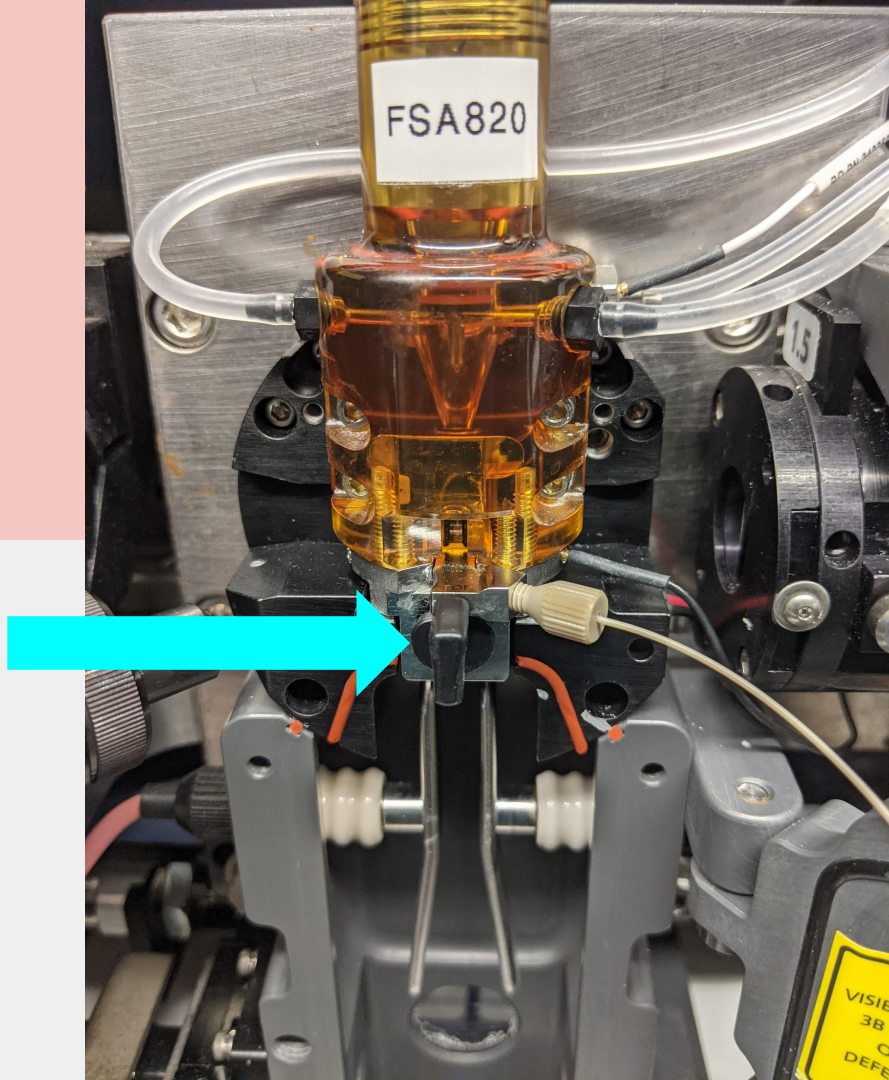

Continue with
the Clean Flow
Cell commands.

Flow Core Staff

Any questions email flowcore@mit.edu